

## ***Moisture determination of animal feed sample***

---

This information is intended as a guideline for development of a moisture analysis method and parameter setting.

Materials vary throughout the world and we will be more than happy to test customers' individual samples.



### **Precisa moisture analysis XM60/120 Halogen**

Description:	powdered animal feed product
Utensils:	aluminium pan
Preparation:	spread thinly on the pan
Sample weight	1.5 g
Temperature program:	standard
Temperature:	105 °C
Stop-Criteria:	Autostop 0.2 % / 60 s
moisture: (average of 4 measurements)	10.15 %
Standard Deviation (+/-):	0.10 %
Average analysis time:	7 min
Reference:	